



PH105 pH Electrode



Installation Instructions

1. Turn meter power "off."
2. Remove retaining collar holding probe on meter.
3. Pull, with a gentle "rocking" motion, to remove old probe.
4. Reverse order to install new probe. Do not touch contacts.
5. Before first use or after extended storage, soak the electrode (with its cap removed) in a pH 4 solution for about 10 minutes.

For further instructions, see instrument User Manual.

This probe is compatible with:

CL200 ExStik Chlorine Meter
PH100 ExStik pH Meter
PH110 ExStik pH Meter
RE300 ExStik ORP Meter

Contact Customer Support

Customer Support Telephone List: <https://support.flir.com/contact>

Calibration, Repair, and Returns: repair@extech.com

Technical Support: <https://support.flir.com>

Copyright © 2022 Teledyne FLIR LLC

All rights reserved including the right of reproduction in whole or in part in any form
www.extech.com